

To whom it my concern,


We wish to inform you, that delivered microwave autoclaves Microjet are equipped with valid software version to allow liquid media sterilization with volume of 500 ml. Delivered manual operating contains incorrect vessel data, which can be use for sterilization.

Please note to mentioned below table with correct range of vessels and mediums to avoid boil over during sterilization process.

We recommended for 500 ml liquid medium to use standard glass bottles with capacity 750 ml or Erlenmayer flask with range capacity from 750 ml up to 800 ml.

TABLE 1. RECOMMENDED STERILIZATION VOLUMES

Container - nominal capacity [ml]	Minimal volume [ml]	Maximum volume [ml]
100	100	100
200	100	150
250	150	200
500	150	400
750	250	500

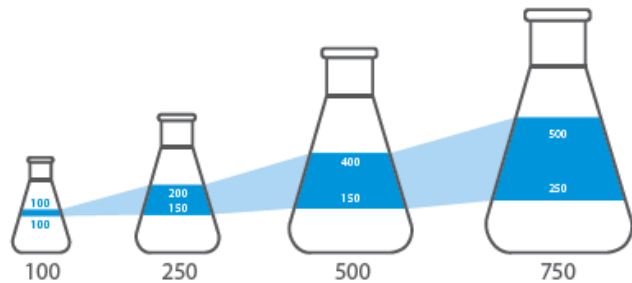

For any further questions please don't hesitate to contact us.

With sincerely

ENBIO team